



**Instructions for the
collection of Tooth Samples**

Kit Contents:

- A 50ml Tooth Container with 25ml of Tooth Transport Solution
- A Preservation Tablet
- Zip lock bag with absorbent towel
- Documentation
- Screw top container

Please follow the instructions listed below:

- 1 When the Tooth has been removed, complete the label on the Tooth Container.
- 2 Dissolve the preservation tablet in the Tooth Transport Solution as follows:



- a) Remove the lid from the Tooth Container. Be careful to place it where it will not be knocked over.



- b) Remove the lid from the preservation tablet container.
- c) Tip the tablet into the Tooth Container and replace the lid.



- 3** The tablet will initially float. Allow the tablet to dissolve - this will take a few minutes.



- 4** As the tablet dissolves, you will see that it sinks to the bottom of the tube.



- 5** Shake the tube well to ensure that the tablet has been evenly distributed through the Tooth Transport Solution.



- 6** Place the tooth in the tooth container.
- Close the lid tightly, ensuring that it is not cross-threaded.



- 7 Place the tube in the zip lock bag with the absorbent towel and place in the screw top container, in the transport box.
- 8 Complete the Documentation and place in the transport box.
- 9 Arrange for the collection as detailed overleaf.

Important things to note

- If you do realise that you've forgotten to add the tablet to the Tube, you can add it after the tooth has been placed in the tube.
- **Do not** put the tablet into anything other than the Tooth Container.
- **Do not** dispose of any of the Tooth Transport Solution.
- **Do not** open the Tooth Container before the tooth has been collected.
- **Do not** add the preservation tablet to the Tooth Container before the tooth is available.
- **Do not** clean the Tooth.

Please contact Future Health on

0800 954 5335

between the hours of

08.00hrs & 20.00hrs

and we will arrange for the courier to collect the sample.
Whilst you are waiting for the courier to collect the sample,
you can place the Tooth Container (which contains the Tooth)
in the refrigerator.

Future Health BIOBANK

10 Faraday Building, Nottingham Science & Technology Park,
University Boulevard, Nottingham NG7 2QP, United Kingdom

Tel: +44 115 967 7707 Fax: +44 115 967 7836

info@fhbb.com

www.fhbb.com